

Command:AFINFO

MS2000 and RM2000 Syntax

Format	AFINFO
Minimum Firmware Version Required	v8.7+

Tiger Syntax

Format	[Addr#]AFINFO
Type	Card-Addressed
Minimum Firmware version Required	v3.26+

This command returns all the values of variables and constants that control Autofocus. It also returns maximum focus value found during last autofocus run, original location and location after frame offset was applied.

Example

```
AFINFO
Best Focus:587
Position Preoffset:  0.0002 mm Afteroffset:  -0.0003 mm
Speed :  1    [AF X]
Travel:0.025000 [AF Y]
Frame Offset:3.500000 [AFC Y]
Hill Offset:70 [AF F]
Contrast:10 [AFC X]
Window Size X:98 Y:98 [AL X Y]
Zero ADJ X:50 Y:90 [AFADJ X Y]
ADC Gain:0 [AFADJ Z]
<LF>
```

Related commands are shown inside square brackets, which you can use to change parameters, where applicable

Sometimes because of limited storage space a lite version of the command may be available. It will only displays Best Focus, Positions preoffset and afteroffset.

```
Best Focus:587
Position Preoffset:  0.0002 mm Afteroffset:  -0.0003 mm
```

[commands](#), [ms2000](#), [autofocus](#)

From:

<https://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<https://www.asiimaging.com/docs/commands/afinfo>

Last update: **2023/08/31 02:00**

