

Command:AFMOVE (AM)

MS2000 and RM2000 Syntax

Shortcut	AM
Format	AFMOVE [X=0 or 1] [Y=0 or 1]
Units	integer code (0 or 1)
Remembered	Using SS Z

Tiger Syntax

Shortcut	AM
Format	[Addr#]AFMOVE [X=0 or 1] [Y=0 or 1]
Units	integer code (0 or 1)
Type	Card-Addressed
Remembered	Using [Addr#]SS Z



At this time this Autofocus 2 feature isn't supported on Tiger controllers.

A mode command that influences subsequent commanded moves. If X=1, then upon completion of a commanded XY move ([MOVE](#) and [MOVREL](#)), for example, MOVE X=123 Y=456 , a multi-axis controller will then automatically initiate an auto-focus.

Y=1 enables the [SAFE_TURRET](#) module. Y=0 disables it.

Response

:A or Error Reply.

Example

```
AM X=1
:A<CR><LF>
```

```
AM X?
:A X=1
```

[commands](#), [autofocus](#), [ms2000](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://asiimaging.com/docs/commands/afmove>

Last update: **2024/03/15 18:18**

