

Command:CDATE (CD)

MS2000 or RM2000 syntax

Shortcut	CD
Format	CDATE
Type	Card-Addressed

Tiger syntax

Shortcut	CD
Format	[Addr#]CDATE
Type	Card-Addressed

This command returns the date and time that the current firmware was compiled.

MS2000 Example

```
CD
Dec 19 2008:16:19:59
```

Tiger Example

```
1CD
Dec 19 2008:16:19:59
```

This example shows that the firmware running was compiled on December 19th year 2008 at 4:19:59 PM.

[commands](#), [tiger](#), [ms2000](#)

From:

<http://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://www.asiimaging.com/docs/commands/cdate>

Last update: **2024/12/02 16:37**

