

Command:HALT (\)

MS2000 or RM2000 syntax

Shortcut	\ (the backslash character)
Format	HALT

Tiger syntax

Shortcut	\ (the backslash character)
Format	HALT
Type	Broadcast or Card-Addressed Command

This command will stop all active motors and other actuators too. If there are no errors, a positive reply of :A will be returned. If the HALT command is given while a commanded move is in motion, the controller will reply with the :N-21 error.

Additional Notes for Tiger cards

It's usually a Broadcast Command but can be used as a Card-Addressed Command as well. When addressed to a specific card, it stops motion on that card only. Note that to use as a Card-Addressed Command the full command HALT must be used instead of the shortcut \, because \ is handled quickly in the command parser.

[commands](#), [tiger](#), [ms2000](#)

From:

<http://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://www.asiimaging.com/docs/commands/halt>

Last update: **2025/04/03 19:16**

