

Command:KI

MS2000 or RM2000 syntax

Shortcut	KI
Format	KI [Axis]=### ...
Units	integer
Remembered	Using SS Z

Tiger syntax

Shortcut	KI
Format	KI [Axis]=### ...
Units	integer
Type	Axis-Specific
Remembered	Using [addr#]SS Z

Sets the servo integral error term constant, the integer value ki. Larger values of ki reduce the time for small errors to be corrected at the finish of a move, but decreases stability if set too large.

MOST USERS DO NOT NEED TO USE THIS FUNCTION!

[commands](#), [tiger](#), [ms2000](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://asiimaging.com/docs/commands/ki>

Last update: **2023/08/31 02:36**

