

Command:LCD

Supported on MS2000 only

Shortcut	LCD
Format	LCD "string"
Units	ASCII Characters

Displays the quoted string on the bottom line of the LCD in place of the version information (DIP SWITCH #2 DOWN).

The quotes are syntactically required. If the quotes are not present, or one is missing, then the command will clear the LCD line. This command is write-only and will not return the contents of the LCD line if queried.

LCD "my super string" causes the last line of the LCD to be updated to: my super string

However, if issued this command: LCD my serial string or LCD ? or just LCD, then the last line of the LCD display will be cleared (with no characters present).

[commands](#), [ms2000](#)

From:

<http://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://www.asiimaging.com/docs/commands/lcd>

Last update: **2025/02/11 19:26**

