

Command:MTIME (MT)

MS2000 or RM2000 syntax

Shortcut	MT
Format	MT [axis] = [time in milliseconds]...
Units	Milliseconds

Tiger syntax

Shortcut	MT
Format	MT [Axis] = [time in milliseconds]...
Units	Milliseconds
Type	Axis-Specific

This command returns the settling time of the last move in milliseconds.

Example

```
MT X
:A 168 0
```

This example shows what is returned after sending the command after a small move.

[commands](#), [tiger](#), [ms2000](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://asiimaging.com/docs/commands/mt>

Last update:

2024/06/13 16:20

