

# Command:SAVESET (SS)

MS2000 or RM2000 syntax

<b>Shortcut</b>	SS
<b>Format</b>	SAVESET [X][Y][Z]

Tiger syntax

<b>Shortcut</b>	SS
<b>Format</b>	[addr#]SAVESET [X][Y][Z]
<b>Type</b>	Card-Addressed

SAVESET allows the user to save current parameters settings to Flash memory.

SAVESET Z, saves settings to flash memory

SAVESET Y, restores previously saved settings after a SAVESET X

SAVESET X, will reload factory defaults upon next power-up

## Reply

Upon the start of execution of this command, the controller will reply with a : .

When the execution is complete, an A will follow the colon.

**Note 1:** During the time interval between the : and the A , no serial or manual moves should be given.

**Note 2:** In MS2000 Versions 6.1u and later (see VERSION command), limit settings (see SETLOW and SETUP) are saved if and only if the SAVEPOS command is issued after the command SAVESET Z.

## Example

MS2000 example

```
SS Z
:A
```

Save the current settings to flash memory.

Tiger example

```
1SS Z
:A
```

Save the Tiger Card at Card Address 1's current settings to flash memory.

[commands](#), [tiger](#), [ms2000](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://asiimaging.com/docs/commands/saveset>

Last update: **2025/03/26 14:33**

