

Command:STOPBITS (SB)

Shortcut	SB
Format	STOPBITS [X?] [X=n]
Units	Integer [1 or 2]
Remembered	Using SS Z

N: Sets the number of stop bits, n, to be used for RS-232 serial communication. The default is one 1 stop bit; the other option is two 2 stop bits. Use the [SAVESET Z](#) command to retain the new stop bit setting after power off.



This command is only available on the MS-2000.

[commands](#), [ms2000](#)

From:

<http://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://www.asiimaging.com/docs/commands/stopbits>

Last update: **2025/05/21 20:05**

