

Command:VERSION (V)

MS2000 or RM2000 syntax

Shortcut	V
Format	VERSION [T]

Tiger syntax

Shortcut	V
Format	[addr#]VERSION
Type	Card-Addressed

This commands returns the current firmware version.

Note: The MS-2000 firmware version no longer contains a character at end, the firmware version after 9.2p is 9.50.

Tiger Example

Firmware version of the card at address 1:

```
1V
:A v3.45
```

MS2000 example

Firmware version of the controller:

```
V
:A Version: USB-9.2m
```

Note: New firmware versions no longer contain a character at the end.

```
V
:A Version: USB-9.52
```

The T parameter outputs the version in Tiger format: *MS-2000 firmware 9.54 or above is required.*

```
V T
:A v9.54
```

This can be useful for software that wants to reuse version parsing and comparison code for both controllers.

[commands](#), [tiger](#), [ms2000](#)

From:

<https://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<https://asiimaging.com/docs/commands/version>

Last update: **2025/03/11 18:05**

