

Speed Truth Firmware Modules

Speed Truth Module

The SPEED_TRUTH firmware module is only used on the MS-2000 controller.

When SPEED_TRUTH is included in the firmware build, the [SPEED](#), [AC](#), and [INFO](#) commands will output the actual values for speed and acceleration rather than the requested values.

Speed Untruth Module

The SPEED_UNTRUTH firmware module is only used on the TG-1000 controller.

This module is not included in any Tiger firmware builds, the default behavior of Tiger is to always report the actual values of speed and acceleration with the SPEED, ACCEL, and INFO commands.

If for some reason you need the old MS-2000 behavior on a Tiger controller, this module exists, but is never used in practice.

[tiger](#), [ms2000](#)

From:

<https://www.asiimaging.com/docs/> - Applied Scientific Instrumentation

Permanent link:

https://www.asiimaging.com/docs/speed_truth_module

Last update: **2025/03/18 19:30**

