

Zeiss Axiovert Observer stage mounting screw locations



Location of three 4 x 12 mm screws that secure stage to Axiovert /AxioObserver. There are two in the front & one in rear. Use 3mm Allen wrench to remove screws & then remove stage.

[zeiss](#), [xystage](#), [axiovert](#)

From:

<https://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

https://www.asiimaging.com/docs/zeiss_axiovert_stage_install

Last update: **2021/09/23 17:15**

