


Warning for Zeiss XY Stage

Installing screws incorrectly Will Damage Stage & Void your Warranty. Follow These Instructions



Use M4x18 mm (M4 x 14mm for AxioImager) Allen Head screws for the rear mounting holes.

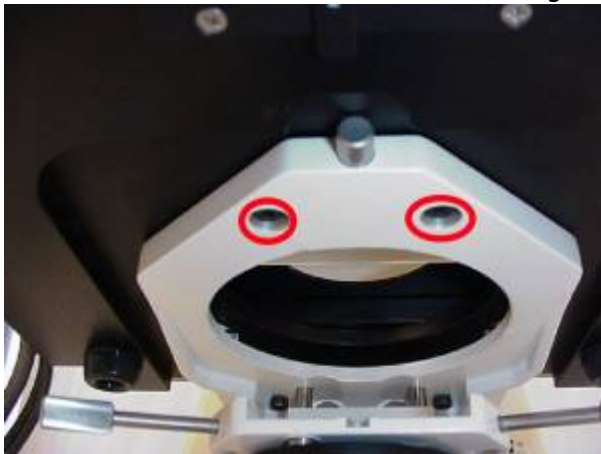

Use M4x10 mm Allen head screws for the front mounting holes.



Before attaching the stage, verify that the mounting screws protrude above the stage carrier no more than 4mm as shown here

If you do not have the correct screws or your stage carrier is different than the one shown here,

please contact ASI before proceeding.



Depending upon the stage carrier you may also need to remove the silver knob shown here by unscrewing it if it interferes with the stage sitting flush / flat on the stage carrier

[xystage](#), [zeiss](#)

From:

<https://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

https://www.asiimaging.com/docs/zeiss_xystage_warning

Last update: **2021/09/23 17:15**

