

## Inserting Pipette into Pipette Holder

1. Disassemble the O-ring compression subassembly from the pipette support subassembly.
2. Insert the pipette through the support subassembly, and through the O-ring.
3. Lightly tighten the outside nut to clamp the pipette without crushing it.
4. Begin screwing the O-ring compression subassembly into the pipette support subassembly. See the two figures below showing the uncompressed O-ring.



O-ring Uncompressed



O-ring Uncompressed

5. Gently tighten the assembly until the O-ring is compressed around the pipette shaft. Be careful not to over-tighten. See the two figures below.



Compressed O-ring



Compressed O-ring