

Command:ZERO (Z)

MS2000 or RM2000 syntax

Shortcut	Z
Format	ZERO

Tiger syntax

Shortcut	Z
Format	ZERO
Type	Broadcast Command

Writes a zero to the position buffer of **all** axes. Allows the user to set current position as the origin. It's a Broadcast Command.

For setting the position of a single axis to zero instead, use the HERE command.

Reply

If there are no errors, a positive response of :A will be returned.

Note

This command has no effect on clocked devices

Example

```
Z  
:A
```

After the reply, the indicators on the LCD should all be zeros.

[commands](#), [tiger](#), [ms2000](#)

From:
<https://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:
<https://www.asiimaging.com/docs/commands/zero>

Last update: **2023/08/31 04:10**

